

Patch for Aphelion Dev and Aphelion Lab version 4.4.0 (build 20083)

This patch fixes the following bugs:

- Images included in the report saved from the Report Generation task is not visible when the report is reloaded in Excel.
- Saving image using the function Save As Image (actual size) from the Window menu does not work when an image had been previously saved on removable volume that is no more available.
- Some Camera configurations based on a IDS uEye camera do not work.

How to patch Aphelion Dev/Lab

If you have not already downloaded the patch file:

1) Download the patch zip file for your Aphelion version available at:

↳ <http://www.adcis.net/en/software-download/>

2) Enter the characters displayed in the image then click the Submit button

After downloading the patch file for Aphelion Dev/Lab 4.4.0 rev 20083:

1) Extract 3 files included in the zip file (Adcis.UI.1.dll, AphelionDev.exe or AphelionLab.exe depending on your Aphelion version, and report_template.xls).

2) Copy these 3 files into the following folder:

↳ For Aphelion Dev 64-bit version: <Aphelion installation folder>\Dev\bin\x64\release

↳ For Aphelion Lab 64-bit version: <Aphelion installation folder>\Lab\bin\x64\release

By default, the <Aphelion installation folder> is:

↳ For 64-bit version: C:\Program Files\ADCIS\Aphelion 4.4.0

These files should replace the files originally installed by the Aphelion Dev/Lab 4.4.0 setup program to correct these issues.